

NCL Method GTA-14

Hep G2 Hepatocarcinoma Homogeneous Apoptosis Assay

Nanotechnology Characterization Laboratory Frederick National Laboratory for Cancer Research Leidos Biomedical Research, Inc. Frederick, MD 21702 (301) 846-6939 ncl@mail.nih.gov

https://ncl.cancer.gov



Method written by:

Stephan T. Stern, Ph.D.

Barry W. Neun, B.S.

Please cite this protocol as:

Stern ST, Neun BW, NCL Method GTA-14: Hep G2 Hepatocarcinoma Homogeneous apoptosis Assay. https://ncl.cancer.gov/resources/assay-cascade-protocols DOI: 10.17917/95VC-ZP26

1. Introduction

This protocol describes the monitoring of nanoparticle treated human hepatocarcinoma cells (HepG2) for apoptosis, as part of the *in vitro* NCL preclinical characterization cascade (1, 2). The protocol utilizes a fluorescent method to measure Caspase-3/7 activation. This is a homogeneous assay that does not require cell isolation, as required for HepG2 Hepatocarcinoma Apoptosis Assay (NCL Method GTA-6).

2. Principles

Caspase-3/7 Homogeneous Fluorometric Assay:

Apoptosis in mammalian cells is initiated by activation of the Caspase family of cysteine proteases. This assay quantifies Caspase-3/7 activation *in vitro* by measuring the cleavage of the non-fluorescent substrate Z-DEVD-R110 (rhodamine 110, bis-(N-CBZ-L-aspartyl-L-glutamyl-L-valyl-L-aspartic acid amide)) to free rhodamine 110, which emits green fluorescence (λ_{max} = 521 nm) (3). This free rhodamine 110 is measured using a microtiter plate reader.

3. Reagents, Materials, Cell Lines, and Equipment

Note: The NCL does not endorse any of the suppliers listed below; their inclusion is for informational purposes only. Equivalent supplies from alternate vendors can be substituted.

- 3.1 Reagents
 - 3.1.1 Acetaminophen (APAP) (Sigma, A7085)
 - 3.1.2 RPMI 1640 (Hyclone, SH30096.01)
 - 3.1.3 L-glutamine (Hyclone, SH30034.01)
 - 3.1.4 Fetal Bovine Serum (Hyclone, SH30070.03)
 - 3.1.5 Nanoparticle
 - 3.1.6 Apo-ONE® Homogeneous Caspase-3/7 Assay Kit (Promega, G7790)
- 3.2 Materials
 - 3.2.1 Costar 96 well flat bottom cell culture plates (Costar, 3598)
- 3.3 Cell Lines
 - 3.3.1 HepG2 (human hepatocarcinoma cells) (ATCC, HB-8065)

3.4 Equipment

- 3.4.1 Plate reader (Safire²-Tecan or equivalent)
- 3.4.2 Plate Shaker

4. Reagent and Control Preparation

- 4.1 Positive control
 - 4.1.1 Acetaminophen (APAP) positive control: Add 8 mg to a total volume of 5 mL RPMI 1640 cell culture media (with 2 mM L-glutamine and 10% FBS) to make a 10 mM solution. Sterile filter using a 0.2 μm filter.
- 4.2 Solutions to make up fresh for each run (can be stored at 4°C for 24 h)
 - 4.2.1 Retrieve Apo-ONE® Assay kit from -20°C.
 - 4.2.2 Thaw the 100X substrate and buffer to room temperature.
 - 4.2.3 Mix by inversion or vortexing. Dilute substrate 1:100 in buffer.
 - 4.2.4 Apo-ONE® Reagent is light sensitive; store protected from light.

5. Experimental Procedure

- 5.1 Cell Preparation (or as recommended by supplier)
 - 5.1.1 Harvest cryopreserved cells from prepared flasks (limit to 20 passages) (Figure 1).
 - 5.1.2 Count cell concentration using a coulter counter or hemocytometer.
 - 5.1.3 Dilute cells to a density of 2.0 x 10⁵ cells/mL in RPMI 1640 cell culture media (2 mM L-glutamine, 10% FBS).
 - 5.1.4 Plate 100 μL/well of diluted cells as per plate format (see Appendix) for each 96 well plate (2 particles/96 well plate).
 - 5.1.5 Incubate plates for 24 h at 5% CO₂, 37°C and 95% humidity (cells should be approximately 80% confluent).
 - 5.1.6 Replace cell culture media with 100 μL media containing test material or positive control. Desired test sample concentration is determined from HepG2 Hepatocarcinoma Cytotoxicity Assay (NCL Method GTA-2). Treat cells for 24 h.

5.2 Caspase Activation Assay

- 5.2.1 Add 100 μ L of Apo-ONE® reagent to each well for a final volume of 200 μ L/well.
- 5.2.2 Cover plate with foil and place on plate shaker (300-500 rpm).
- 5.2.3 Incubate for 1 h at room temperature.
- 5.2.4 Read fluorescence on plate reader with an excitation of 499 nm and an emission of 521 nm.

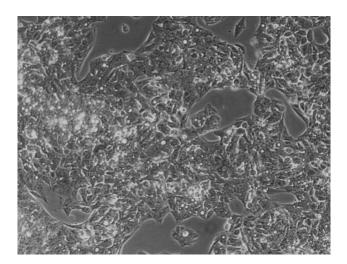


Figure 1. Example of HepG2 Cell Culture Appearance.

Image was taken with a phase contrast microscope at 225X magnification. Human hepatocarcinoma cells (HepG2) are approximately 80% confluent at this stage.

6. Calculations

 $Total\ Caspase\ Activity = \frac{(sample\ fluorescence\ -\ sample\ cell\ free\ blank\ fluorescence)}{mean\ of\ (media\ fluorescence\ -\ media\ blank\ flourescence)} \ *100$

Mean, SD and %CV should be calculated for each positive control and sample.

7. Acceptance Criteria

- 7.1 If a concentration responsive increase or decrease in sample cell-free blank fluorescence is observed, then the nanoparticle is interfering with the assay fluorescence and the non-homogeneous apoptosis assay; NCL Method GTA-6 Hepatocarcinoma Apoptosis Assay should be used instead.
- 7.2 The fold increase in caspase activity at 24 h for the APAP positive control versus media negative control should be at least 3 (300% of negative control).
- 7.3 The positive and sample replicate coefficient of variations should be within 50%.
- 7.4 The assay is acceptable if condition 7.1 and 7.2 are met. Otherwise, the assay should be repeated until acceptance criteria are met.

8. References

- 1. ISO 10993-5 Biological evaluation of medical devices: Part 5 Tests for *in vitro* cytotoxicity.
- 2. F1903 98 Standard Practice for Testing for Biological Responses to Particles *in vitro*.
- 3. Promega Technical Bulletin #295: Apo-ONE® Homogeneous Caspase-3/7 Assay.

9. Abbreviations

APAP acetaminophen

CV coefficient of variation

FBS fetal bovine serum

rpm revolutions per minute

RPMI Rosewell Park Memorial Institute

SD standard deviation

Z-DEVD-R110 Rhodamine 110, bis-(N-CBZ-L-aspartyl-L-glutamyl-L-valyl-L-

aspartic acid amide

10. Appendix

Example of a 96-well plate template.

	1	2	3	4	5	6	7	8	9	10	11	12
	Cells	Cells	Cells	Cells	Cells	Cells	Cells	Cells	Cells	Cells	Cells	Cells
A		Samp. 1	Samp. 1	Samp. 1	Samp. 1	Samp. 1	Samp. 1	Samp. 1	Samp. 1	Samp. 1	Samp. 1	APAP,
	Media	Dilution 10	Dilution 9	Dilution 8	Dilution 7	Dilution 6	Dilution 5	Dilution 4	Dilution 3	Dilution 2	Dilution 1	10 mM
	Cells	Cells	Cells	Cells	Cells	Cells	Cells	Cells	Cells	Cells	Cells	Cells
В		Samp. 1	Samp. 1	Samp. 1	Samp. 1	Samp. 1	Samp. 1	Samp. 1	Samp. 1	Samp. 1	Samp. 1	APAP,
	Media	Dilution 10	Dilution 9	Dilution 8	Dilution 7	Dilution 6	Dilution 5	Dilution 4	Dilution 3	Dilution 2	Dilution 1	10 mM
	Cells	Cells	Cells	Cells	Cells	Cells	Cells	Cells	Cells	Cells	Cells	Cells
C		Samp. 1	Samp. 1	Samp. 1	Samp. 1	Samp. 1	Samp. 1	Samp. 1	Samp. 1	Samp. 1	Samp. 1	APAP,
	Media	Dilution 10	Dilution 9	Dilution 8	Dilution 7	Dilution 6	Dilution 5	Dilution 4	Dilution 3	Dilution 2	Dilution 1	10 mM
	Media	Media	Media	Media	Media	Media	Media	Media	Media	Media	Media	Media
D		Samp. 1	Samp. 1	Samp. 1	Samp. 1	Samp. 1	Samp. 1	Samp. 1	Samp. 1	Samp. 1	Samp. 1	APAP,
	Media	Dilution 10	Dilution 9	Dilution 8	Dilution 7	Dilution 6	Dilution 5	Dilution 4	Dilution 3	Dilution 2	Dilution 1	10 mM
	Media	Media	Media	Media	Media	Media	Media	Media	Media	Media	Media	Media
E		Samp. 1	Samp. 1	Samp. 1	Samp. 1	Samp. 1	Samp. 1	Samp. 1	Samp. 1	Samp. 1	Samp. 1	APAP,
	Media	Dilution 10	Dilution 9	Dilution 8	Dilution 7	Dilution 6	Dilution 5	Dilution 4	Dilution 3	Dilution 2	Dilution 1	10 mM
	Cells	Cells	Cells	Cells	Cells	Cells	Cells	Cells	Cells	Cells	Cells	Cells
F		Samp. 2	Samp. 2	Samp. 2	Samp. 2	Samp. 2	Samp. 2	Samp. 2	Samp. 2	Samp. 2	Samp. 2	APAP,
	Media	Dilution 10	Dilution 9	Dilution 8	Dilution 7	Dilution 6	Dilution 5	Dilution 4	Dilution 3	Dilution 2	Dilution 1	10 mM
	Cells	Cells	Cells	Cells	Cells	Cells	Cells	Cells	Cells	Cells	Cells	Cells
G		Samp. 2	Samp. 2	Samp. 2	Samp. 2	Samp. 2	Samp. 2	Samp. 2	Samp. 2	Samp. 2	Samp. 2	APAP,
	Media	Dilution 10	Dilution 9	Dilution 8	Dilution 7	Dilution 6	Dilution 5	Dilution 4	Dilution 3	Dilution 2	Dilution 1	10 mM
	Cells	Cells	Cells	Cells	Cells	Cells	Cells	Cells	Cells	Cells	Cells	Cells
Н		Samp. 2	Samp. 2	Samp. 2	Samp. 2	Samp. 2	Samp. 2	Samp. 2	Samp. 2	Samp. 2	Samp. 2	APAP,
	Media	Dilution 10	Dilution 9	Dilution 8	Dilution 7	Dilution 6	Dilution 5	Dilution 4	Dilution 3	Dilution 2	Dilution 1	10 mM